WALL MOUNTED – WANDMONTAGE – VÆGMONTAGE

UK – System 50 + 63 + 75 WM

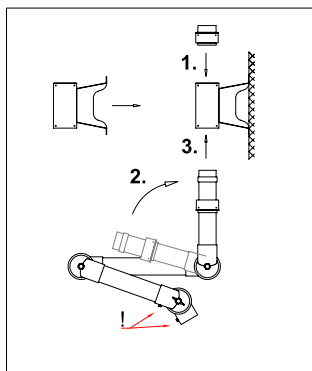
Fig. 1

1. Fasten the wall bracket solidly to the wall or column. When mounting the wall bracket be aware of the considerable strain the operation of the arm will transmit to the bracket.
2. Place the square-guide in the top of the bracket (1) and fasten it to the wall bracket with the included screws.
3. Considering the internal spring, it is important to turn the socket pipe in the direction of the arrow (2) and that the fingerscrews (!) are faced forward.
4. Place the socket pipe in the wall bracket(3) and fasten the square-guide with the included screws.
5. The extraction arm can be turned 190° in the wall bracket.

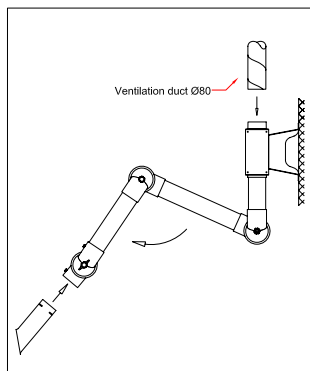
Fig. 2

1. Connect the main ventilation duct (80 mm dia.) directly to the top of the wall bracket.
2. Unfold the extraction arm in the direction of the arrow.
3. Place the connection pipe of the accessory in the joint. Turn the slit on the connection pipe towards the small fingerscrew on the joint. When screwed in the fingerscrew will hold the accessory in place in the joint.
4. Finally make sure the damper is open.
5. The extraction arm is now ready for use.

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See MAINTENANCE AND CLEANING
at www.alsident.com under "Technical Information"

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UK – System 50 + 63 + 75 CM

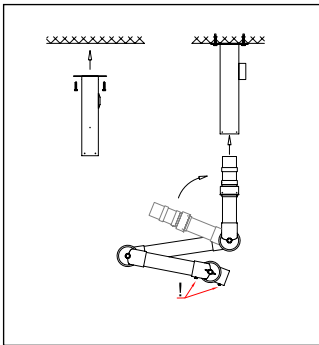
Fig. 1

1. Mount the column under the ceiling.
2. The operation of the extraction arm will transmit a considerable strain to the column. The column must therefore be mounted on a solid ceiling.
3. Considering the internal spring, it is important to turn the socket pipe in the direction of the arrow and that the fingerscrews (!) are faced forward.
4. Place the socket pipe in the ceiling column and fasten the square-guide with the included screws.
5. The extraction arm can now be turned 360° in the column.

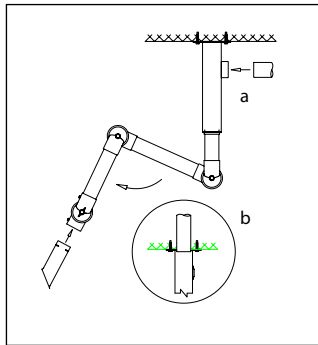
Fig. 2

1. For side connection (Fig. 2a) the main ventilation duct (80mm dia.) can be connected directly to the outlet on the side of the column.
For top connection (Fig. 2b) the main ventilation duct is put through the ceiling and directly into the column. Make sure to tighten the gap between the ceiling and duct with a filler.
2. Unfold the extraction arm in the direction of the arrow.
3. Place the connection pipe of the accessory in the joint. Turn the small slit in the connection pipe towards the fingerscrew. When screwed in the fingerscrew will hold the accessory in place in the joint.
4. Finally make sure the damper is open.
5. The extraction arm is now ready for use.

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All rights reserved
Alsident® System A/S