
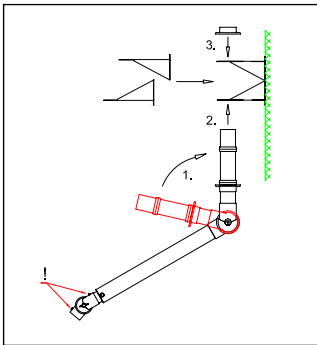
# WALL MOUNTING – WANDMONTAGE – VÆGMONTAGE

## UK – Teleskop – WM

**Fig. 1**


1. The wall bracket is in two parts. Mount the bracket solidly on the wall or column and close together. When mounting the wall bracket be aware of the considerable strain the operation of the arm will transmit to the wall bracket.
2. Considering the internal spring, it is important to turn the socket pipe in the direction of the arrow and that the fingerscrews (!) are faced forward.
3. Place the socket pipe in the bracket and fasten the flange to the bracket with the screws.
4. Place the upper flange on the socket pipe and fasten it to the bracket with the included screws.
5. The two flanges are not fastened to the socket pipe. Therefore the extraction arm can be turned 190° in the wall bracket.

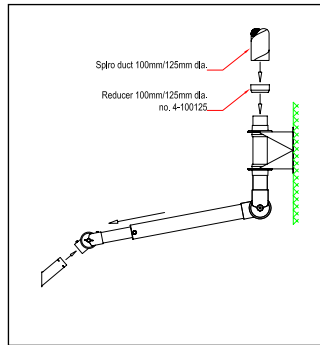**Fig. 2**


1. Loosen the fingerscrew located at the pipe reduction and pull the extraction arm to the length. Tighten the fingerscrew slightly again.
2. Connect the extraction arm to the main ventilation duct with the Alsident® reducer. In order to turn the extraction arm, it is important not to fix the reducer to the socket pipe.
3. Place the connection pipe of the accessory in the joint. Turn the slit on the connection pipe towards the small fingerscrew. When screwed in the fingerscrew will hold the accessory in place in the joint.
4. Finally make sure the damper is open.
5. The extraction arm is now ready for use.

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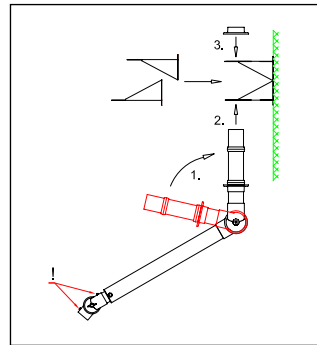
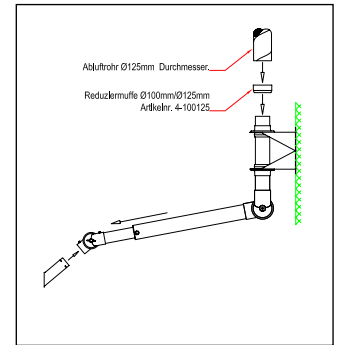
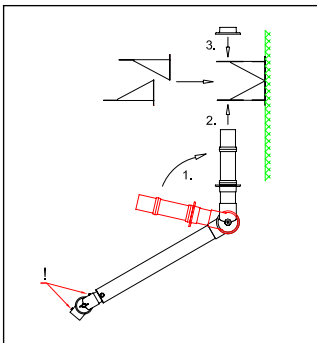
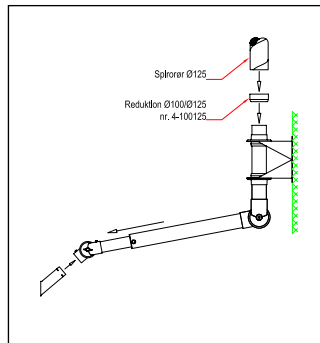
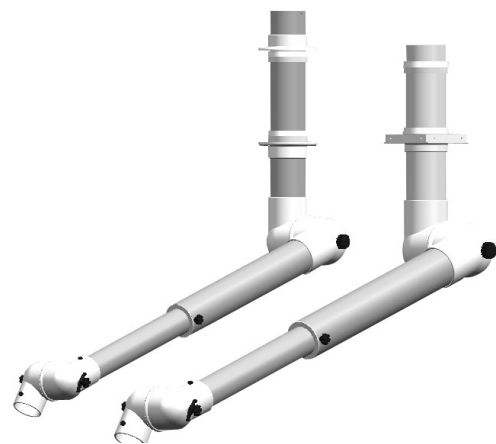
## UK – Teleskop – CM

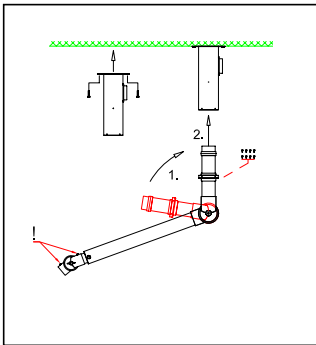**Fig. 1**


1. The operation of the extraction arm will transmit a considerable strain to the column. The column must therefore be mounted on a solid ceiling.
2. Considering the internal spring, it is important to turn the socket pipe in the direction of the arrow and that the fingerscrews (!) are faced forward.
3. Place the socket pipe in the column. Turn the guide on the socket pipe until it falls into place in the column and fasten with the included screws.
4. The extraction arm can now be turned 360° in the column.

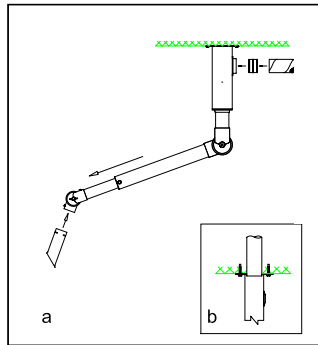**Fig. 2**


1. For side connection (fig. 2a) the main ventilation duct (125 mm dia.) can be connected to the outlet on the side of the column with a nipple – or directly to the outlet.  
For top connection the main ventilation duct is put through the ceiling and directly into the column. Make sure to seal the gap between the ceiling and duct with a filler.
2. Loosen the fingerscrew located at the pipe reduction and pull the extraction arm to the length. Tighten the fingerscrew slightly again.
3. Place the connection pipe of the accessory in the joint. Turn the slit on the connection pipe towards the small fingerscrew. When screwed in the fingerscrew will hold the accessory in place in the joint.
4. Finally make sure the damper is open.
5. The extraction arm is now ready for use.

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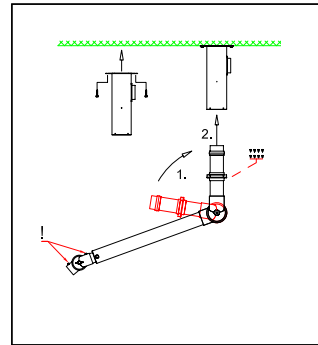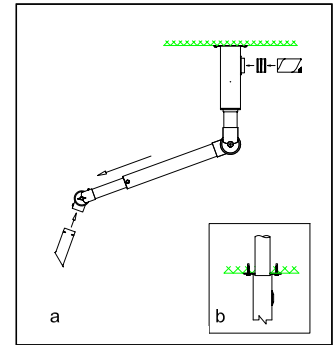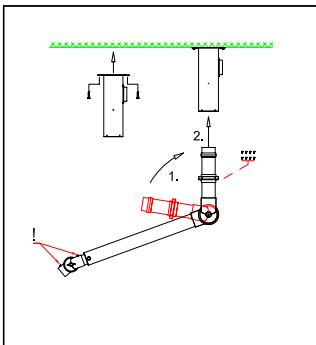
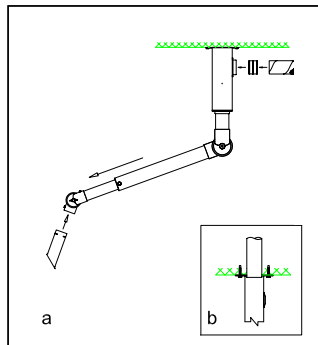
See MAINTENANCE AND CLEANING  
at [www.alsident.com](http://www.alsident.com) under "Technical Information"

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All rights reserved